


UVL-4F Longwave UV Lamp

User's Manual

Important: Please read the instructions carefully prior to using the lamp in order to avoid malfunction or damage.

Operation

Easily switch between the longwave UV and the flashlight by using the selector located on the side of the lamp.

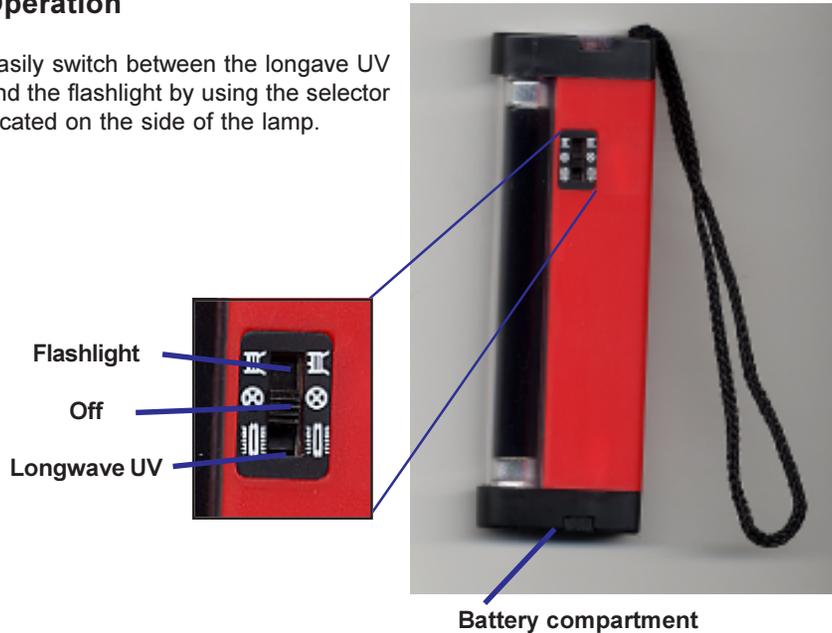

Insert 4 AA batteries (UM-3) not included with the lamp. To insert batteries, slide the battery cover out and flip upwards. Position batteries as shown on the inside of the battery compartment. Slide back the battery cover.

Keep the lamp away from extreme temperature changes, excessive humidity, moisture, rain or other harsh environments. Do not mix old and new batteries together. It is not recommended to use the fluorescent lamp function when the energy level is low, it will blacken the fluorescent tube. Remove the batteries if the lamp is not to be used for a long period.


Replacement Parts
Replacement Tube, 4W longwave BLB

Part Number
34-0010-01



Replacing the Longwave Tube or Bulb

1. Remove the bulb cover. The cover is tightly screwed into the housing. To remove, unscrew in a one-quarter counter-clockwise turn. Lift off the bulb cover and the black end cap.
2. Remove the flashlight bulb as shown in the diagram. Replace with a 4.8V bulb.
3. To remove the longwave tube, slide off the clear tube cover.
4. Turn the tube slightly and it will slide out. Replace with a 4watt longwave tube. Replace the clear cover.
5. Reconnect the end cap. Slide bulb cover back into position, aligning the notches. Turn cover clockwise to close.

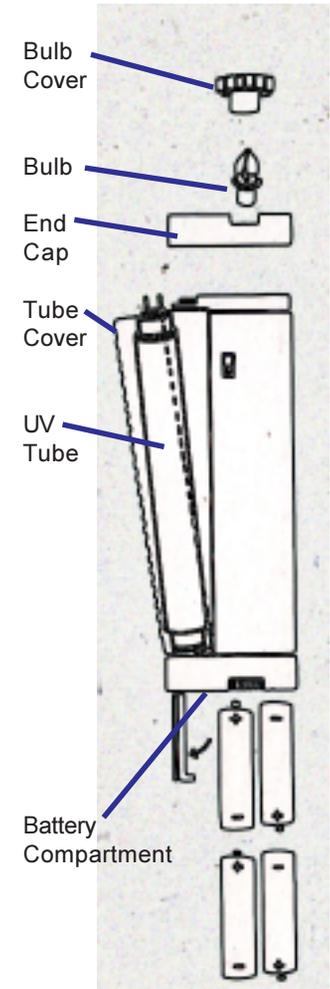

Replacing the Batteries

1. Slide the battery cover out and lift it up.
2. Remove the old batteries and replace with four new "AA" size "UM3" batteries, noting the positive and negative indication mark for proper insertion.
3. Slide the battery cover back into position.



UVP, Inc. 2066 W. 11th Street, Upland, CA 91786 Tel: (800) 452-6788 or (909) 946-3197 Fax: (909) 946-3597
Ultra-Violet Products Ltd. Unit 1, Trinity Hall Farm Estate, Nuffield Road, Cambridge CB4 1TG UK Tel: +44(0)1223-420022 Fax: +44(0)1223-420561
Web Site: www.uvp.com